

Technical Note SBT Resolver™ DRB1 ST-PD5.2-8, Dec-2015

PRODUCT CODE

ST-PD5.2-8(50)

ST-PD5.2-8(20)


UPDATE DESCRIPTION

The new ST-PD5.2-8 assay now contains the DNA polymerase within the PCR mix. This new single-tube formation allows for improved useability and increased assay robustness with PCR uniformity and strong signal intensity.

IMPACT OF UPDATE

There is no longer a DNA-Polymerase vial supplied within the kit. Therefore it is recommended that users carefully check the IFU as the PCR dispense volumes have changed. As a consequence of this, the total volume in the HLA-DRB1 PCR MIX vial has also increased to accommodate the altered dispense volume.

For a copy of the ST-PD5.2-8 IFU, ask your distributor or email us at support@conexio-genomics.com.

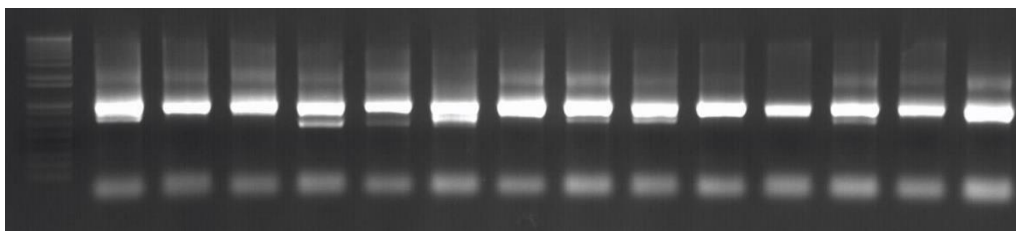

DRB1 ST-PD5.2-8 assay for multiple samples demonstrating uniformity of amplification

ANALYSIS IN ASSIGN™

Analysis for the ST-PD5.2-8 assay in Assign™ remains the same as for the LG-PD5.2-7 assay and uses the same 527_DRB1.xml reference file.

It is recommended that users name EPGs with '8DRB1' to differentiate the assay historically, in which case the alias '8DRB1' should be added to the 527_DRB1.xml reference.



CONEXIO

Pioneering since 2002

The new **HLA-DRB1 ST-PD5.2-8** assay, within the SBT Resolver™ suite, is for **Research Use Only**. Not for use in diagnostic procedures.
No claim or representation is intended to provide information for the diagnosis, prevention or treatment of a disease.

If you have any questions in regard to our new product, please contact your local Olerup representative or contact Conexio Genomics at support@conexio-genomics.com



Conexio Genomics Pty Ltd
2/49 Buckingham Dr
Wangara 6065
Western Australia
Australia

Distributed by



www.olerup.com

Designed, developed and manufactured by Conexio

conexio-genomics.com

TEC257

Version 1.0