I N S T R U C T I O N M A N U A L

***OLERUP* SSP HORIZONTAL
GREEN
SYSTEM**

Model:

Gel System 96

C.B.S. ☆ SCIENTIFIC COMPANY, INC.

P.O. Box 856, Del Mar, CA 92014
619-755-4959 or Toll-free 800-243-4959

TABLE OF CONTENTS

	Page
Important User Information	3-4
Section 1 General Information	
1.1 Introduction	5
1.2 Specifications	5
1.3 Safety	5
Section 2 Description of parts	
2.1 Unpacking	6
2.2 Components/Assembly	6
Section 3 Instructions for Use	
3.1 Gel Unit Preparation and Gel Casting	7-8
3.2 Running the Gel	8
3.3 Removing the Gel	8
3.4 Maximum Well/Comb volumes	8
Section 4 Running Conditions	
4.1 Recommended Power	9
4.2 Recommended Buffers	9
4.3 References	9
Section 5 Maintenance of Equipment	
5.1 Care and Handling	10
5.2 Maintenance	10
Section 6 Equipment and Accessories	11
Addresses	12

IMPORTANT USER INFORMATION

This Instruction Manual will explain how to use this product safely and effectively. Please read and carefully follow the instruction manual in its entirety.

The triangle/exclamation mark symbol alerts the user of the product to important operational, maintenance, and/or warranty requirements.

The triangle/lightning bolt symbol alerts the user of the product to potentially hazardous electrical exposure.

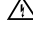
Failure to adhere to the instructions could result in personal and/or laboratory hazards, as well as invalidate any warranty. Always turn off the DC power source prior to disconnecting power cords from the product. Disconnect power cords from the power source first, and then from the product. For maximum safety, always operate this system in an isolated, low traffic area, not accessible to unauthorized personnel. Never operate damaged or leaking equipment.

WARRANTY AND LIABILITY

This product was produced utilizing the highest practical standards of materials, workmanship, and design. C.B.S. Scientific warrants that the product has been tested and will meet or exceed published specifications. This warranty is valid only if the product has been operated and maintained according to the instructions provided.

C.B.S. Scientific warrants this product to be free from defects in materials and workmanship under normal service for one year from date of shipment. If the product proves defective during this period, C.B.S. Scientific will repair or replace it at our option, free of charge, if returned to us postage prepaid. This warranty does not cover: damage in transit, damage caused by carelessness, misuse or neglect, normal wear through frequent use, damage caused by solvent corrosion, damage caused by improper handling or user alteration, nor unsatisfactory performance as a result of conditions beyond our control. C.B.S. Scientific shall in no event be liable for incidental nor consequential damages, including without limitation, lost profits, loss of income, loss of business opportunities, loss of use and other related damages, however caused, nor any damage arising from the incorrect use of the product.

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General Information

1.1 Introduction

The green Horizontal unit, Gel System 96 (16 x 17.5 cm) effectively allows rapid separation and identification of various fragments from DNA, RNA, PCR™ products and synthetic oligonucleotides. When used in combination with two or more sample combs, designed for multi-channel pipette loading, multiple screening of large sample numbers can be facilitated easily.

1.2 Specifications

Constructions:

Buffer chamber	Acrylic
Safety Cover	Acrylic
Electrodes	Platinum wire .012" diameter
Power cords	FR Urethane rated 7500VDC, 200mA, 65°C
Combs	Teflon coated aluminum
UVT Gel Tray	UV transparent acrylic
Safety Certification	EN61010-1-1993 (IEC1010-1)

	Gel System 96
Shipping weight	5 lbs
Overall Size (l)x(w)x(h) cm	28 cm x 21 cm x 13 cm
Gel Size cm	16 cm x 17.5 cm
Recommended buffer volume: with gel tray & gel	750 ml
Distance between electrodes	25 cm
Voltage limit	250VDC

1.3 Safety

Power to the Gel System 96 unit is to be supplied by an external DC voltage power supply that must be ground isolated, so that the DC voltage output floats with respect to ground. For any power supply used, the maximum specified operating parameters for the units are:

Maximum Limits

250 VDC voltage
40 watts power
40 mA current
55°C ambient temperature

Current to the unit, provided from the external power supply, must enter the unit through the lid assembly, providing a safety interlock to the user. Current to the unit is broken when the lid is removed. **Do not attempt to use the unit without the safety lid, and always turn the power supply off before removing the lid, or when working with the unit in any way. Follow safety precautions specified by the power supply manufacturer.**

SECTION 2


Description of Parts

2.1 Unpacking

Please verify that your unit comes complete with the following components:

- Green Horizontal electrophoresis unit
- Safety cover with attached DC power leads
- UV Transparent slotted gel tray
- 6 Aluminum Taper Tooth combs

2.2 Components/Assembly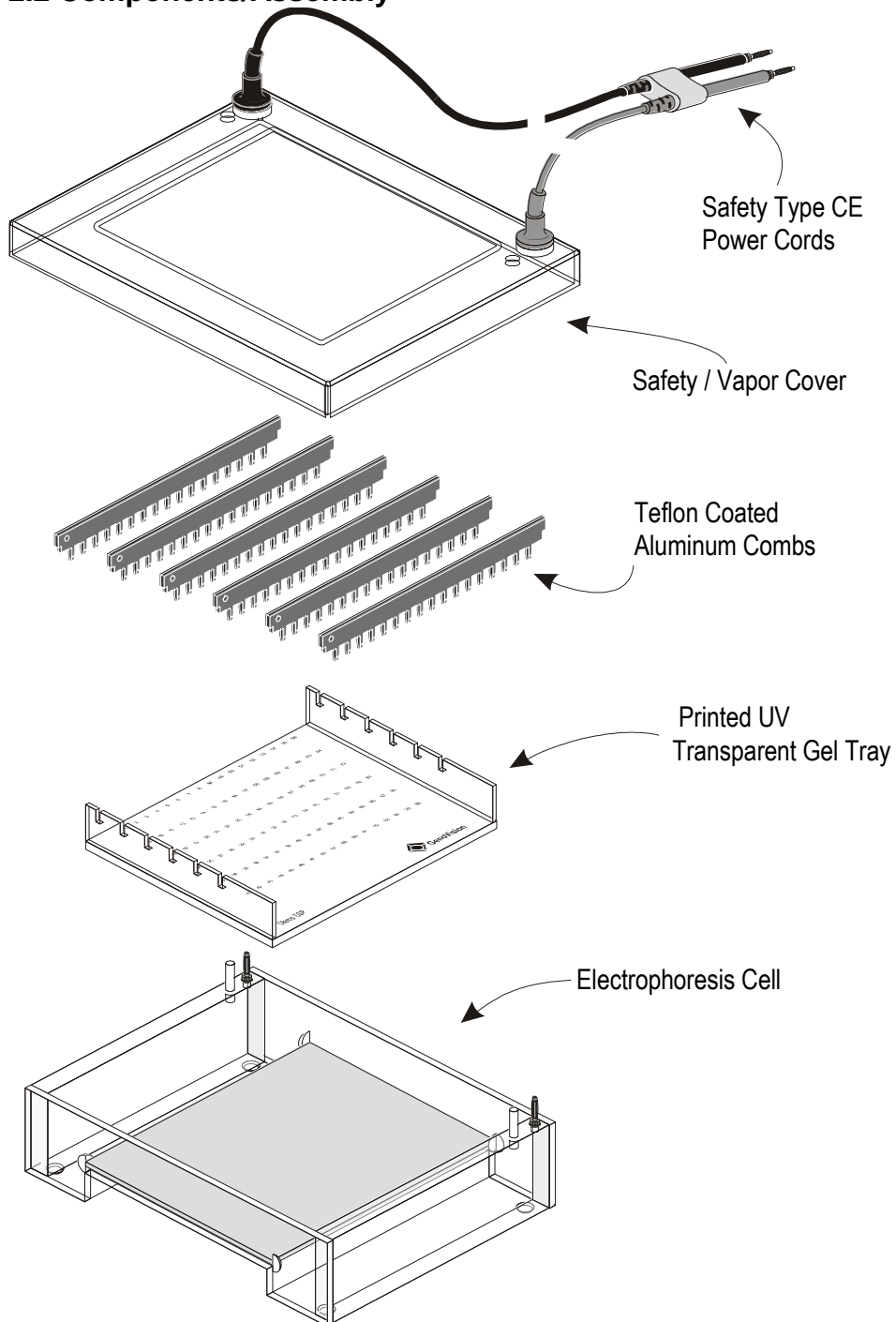

SECTION 3

Instructions for Use

3.1 Gel Unit Preparation and Gel Casting


Place gel unit on a level work surface in an authorized work area. Place the combs in the gel tray. **See Table 1 for comb volumes.**


A. Gel Casting using the Single or Multiple Gel Caster

1. Place the gel casting unit on a lab bench in an approved work area. Level the unit using the white delrin leveling screws and bubble level.

Gel Casting Using Single "Flexcaster"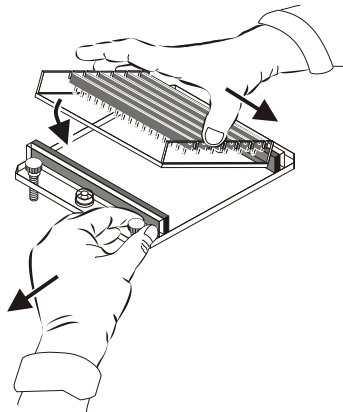

2. Insert the gel tray by pressing one end against the rubber sponge gasket. Quickly lower the other end onto the gel caster and release against other gasket. (**See drawing**). Repeat, if using a multiple gel caster.

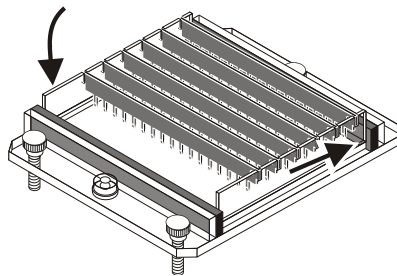3. Insert combs into gel tray slots.

4. Over a Bunsen burner or in a microwave, boil agarose/buffer solution. Recommended: 2% (w/v) agarose in 0.5 x TBE buffer, 100 ml of gel solution per gel tray. Allow to cool to 55-60°C before pouring onto gel tray. Preferably stain the gel prior to casting with ethidium bromide (10mg/ml, 5 µl per 100 ml gel solution. For maximal ease of handling use our ethidium bromide dropper bottles, 1 drop of ethidium bromide solution per 50 ml of gel, Prod. No. 103.301-10).

5. Allow to cool to room temperature. Remove combs and transfer to gel electrophoresis unit.

3.2 Running the gel

1. Add enough buffer to fill both reservoirs and overflow the surface of the gel to a depth of 2-3mm. Gently remove comb(s). Flush out any air bubbles in the wells. Load the samples into the sample wells.

NOTE: DO NOT FORGET TO LOAD DNA SIZE STANDARD.
(100 base pair ladder, Prod. No. 103.201-100.)

2. Align safety cover over the unit and carefully attach, so not to disturb samples.

3. Connect the leads to the power supply, matching the color-coded red to red (the anode, +) and black to black (the cathode, -).

See Section 4.1 for recommended power conditions.

Begin separation by electrophoresis.

3.3 Removing the Gel

1. Turn the power supply off and disconnect the leads from the power supply.

2. Remove the safety cover from the unit, by placing thumbs on white posts next to red & black connectors, then pushing down while pulling up with fingers under lid.

DO NOT pull on power cords.

3. Gently lift the gel tray from the unit. **Always wear gloves, eye protection and protective clothing** if buffer and/or gel contains Ethidium Bromide. Ethidium Bromide is a powerful mutagen, gloves, eye protection and protective clothing should always be worn when handling the gel or buffer solutions. See Material Data Safety Sheets.

4. View separated fragments under UV light, using proper protection for eyes and skin (see manufacturer's instructions).

3.4 Maximum Well/Comb Volumes

Tooth Width (mm)	Tapered Tooth Thickness @ tip	Maximum Volume per 1mm Sample Height
3 mm	0.8 mm	2.4 μ l

Rec. Volume of Sample = 5-10 μ l

SECTION 4 Running Conditions

4.1 Recommended Power

As a rule, optimal resolution of larger molecules is achieved during longer runs at lower voltages, whereas smaller molecules require shorter runs at higher voltages. Applied voltage gradients can therefore be anywhere in the range of 1-10 VDC/cm of gel. Using the standard buffer systems listed below, most runs will use 5 VDC/cm as a general rule. Recommended: 8 VDC/cm.

The usual run time will vary for the voltage chosen but should range from 15 to 60 minutes. Recommended: 20 to 30 minutes. Nucleic acid migration is monitored by the progress of marker dyes. Constant power is not a necessity, but it produces uniform heat throughout the run, therefore minimizing band diffusion. Be sure the polarity is correct i.e. that the DNA is loaded near the cathode (black electrode) to run toward the anode (red terminal).

Agarose gels may be stored for several days at 4°C wrapped in plastic wrap. Seakem Agarose (FMC) is used (normally) for preparative and analytical gels. Other types of agarose can be used for special purposes.

4.2 Recommended Buffers*

Type*	Concentrated Stock/liter	Final Concentration
TAE (Tris-acetate)	50X - 242 gm Tris base 57.1ml glacial acetic acid 100 ml 0.5M EDTA (pH 8.0)	1X - 0.04M Tris-acetate 0.001M EDTA
TBE (Tris-borate)	5X - 54 gm Tris base 27.5 gm boric acid 20 ml 0.5M EDTA (pH 8.0)	0.5X - 0.045M Tris-borate 0.001M EDTA

Ethidium Bromide Staining

EtBr can be premixed with buffers and agarose for use during electrophoresis. Add to agarose only after temperature has fallen below 55°C. Gels can also be stained after electrophoresis in a soaking tray. Use EtBr at a final concentration of 0.1µg/ml from a 1mg/ml stock solution. Ethidium Bromide is a powerful mutagen. ALWAYS wear gloves, eye protection and protective clothing. Dispose of solutions in accordance with the safety regulations of your institution.

4.3 References

1. Hames, B.D., Rickwood, D. (ed.) (1990). *Gel Electrophoresis of Nucleic Acids. A Practical Approach*. 2nd edn. IRL Press, Oxford. Ch. 2.
2. Sambrook, J., Fritsch, E.F., Maniatis, T. (1989). *Molecular Cloning. A Laboratory Manual*. 2nd edn. Cold Spring Harbor Laboratory Press, Cold Spring Harbor, New York. Ch 6.
3. Ausubel, F.M., Brent, R., Kingston, R.E., Moore, D.D., Seidman, J.G., Smith, J.A., Struhl, K. (ed) (1993). *Current Protocols in Molecular Biology*. Vol. 1, Greene Publishing Associates, Inc. and John Wiley & Sons, Inc., Ch. 2.

SECTION 5 Maintenance of Equipment

5.1 Care and Handling

The plastic components of the Midi-Horizontal units are fabricated from acrylic and polycarbonate. Electrodes and connectors are made from pure platinum, stainless steel, and chrome plated brass. As with any laboratory instrument, adequate care ensures consistent and reliable performance.

After each use, rinse buffer chamber, gel tray and combs with de-ionized water. Wipe dry with a soft cloth or paper towel, or allow to air dry. Whenever necessary, all components may be washed gently with water and a non-abrasive detergent, and rinsed and dried as above. *Never* use abrasive cleaners, glass cleaning sprays or scouring pads to clean the components, as these will damage the unit and components.

Additional precautions:

- Do not autoclave or dry-heat sterilize the apparatus or components.
- Do not expose the apparatus or components to phenol, acetone, benzene, halogenated hydrocarbon solvents or undiluted alcohol's.
- Avoid prolonged exposure of the apparatus or components to UV light.
- Do NOT treat with diethylpyrocarbonate (DEPC)-treated water for extended periods at 37°C. A brief rinse with DEPC-water is sufficient after a thorough wash, followed by a quick rinse in 70% ethanol.

5.2 Maintenance

The following inspection and maintenance procedures will help maintain the safety and reliable performance of the Green Horizontal systems. Replacement parts can be ordered by call 1-858-755-4959 or by contacting your local distributor.

- Banana plugs and power cords should be inspected regularly. If the banana plugs become loose or do not feel friction tight replace the plugs or power cords.
- Should power cord assemblies (connectors, wire or shrouds) show any signs of wear or damage (e.g. cracks, nicks, abrasions, or melted insulation), replace them immediately.
- The platinum wire is secured to the banana jack by compression between a stainless washer and the jack nut. The nut/washer interface should be tight and free of corrosion.

SECTION 6

Olerup SSP Horizontal Green System and Accessories

Product No.	Description
103.101-01	Gel System 96. Includes: basic gel unit with NO tapered gates, safety cover with attached power leads, UVT Precision / silk-screened gel tray, 6 each combs (aluminum teflon coated, double tapered)

Gel Trays

Product No.	Description
103.101-11	UVT precision/ silk-screened gel tray

Combs

Product No.	Description
103.101-21	17 wells, double tapered, aluminum teflon coated

Gel Casters

Product No.	Description
103.101-31	Gel caster for one gel. One gel tray included.
103.101-33	Gel caster for three gels. Three gel trays included.

Related products

Product No.	Description
103.301-10	Ethidium bromide dropper bottle. 10 ml. 1 drop per 50 ml of gel solution.
103.302-05	GelRed Dropper Bottle 5 ml, 4 drops for every 100 to 120 ml of gel solution.
103.201-100	DNA size marker. 100 base pair ladder, 10 fragments 100 to 1000 base pairs. 100 applications
103.203-100	DNA Size Marker for short gel runs

ADDRESSES:

Manufacturer:

Olerup SSP AB, Franzengatan 5, SE-112 51 Stockholm, Sweden.

Tel: +46-8-717 88 27

Fax: +46-8-717 88 18

E-mail: info-ssp@olerup.com

Web page: <http://www.olerup-ssp.com>

Distributed by:

Olerup GmbH, Löwengasse 47 / 6, AT-1030 Vienna, Austria.

Tel: +43-1-710 15 00

Fax: +43-1-710 15 00 10

E-mail: support-at@olerup.com

Web page: <http://www.olerup.com>

Olerup Inc., 901 S. Bolmar St., Suite R, West Chester, PA 19382

Tel: 1-877-OLERUP1

Fax: 610-344-7989

E-mail: info.us@olerup.com

Web page: <http://www.olerup.com>

For information on *Olerup* SSP distributors worldwide, contact **Olerup GmbH**.